IDEXX SNAP* MRL Beta-Lactam Test

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Precautions: Test milk within 3 days of collection; thoroughly mix milk before testing. Do not use abnormal-looking milk or expired tests; do not mix components from different kits. If the control spot fails to develop color, retest the sample.

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

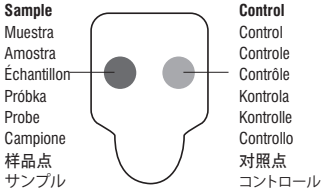
OR
O
OU
OU
LUB
ODER
OPPURE
或
または

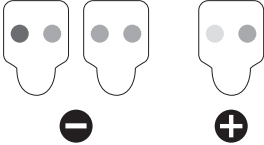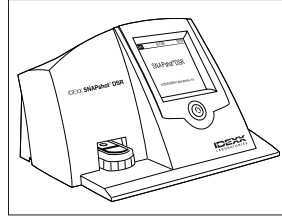IDEXX SNAPshot* DSR Reader

It is recommended that positive samples be further analyzed.

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Technical assistance • idexx.com/dairy

US/Canada/Latinoamerica	+1 800 321 0207	日本	+021 6127 9528
EU	+00800 329 43399	Within AU	+1300 44 33 99
中国	+010 84423133	Within NZ	+0800 102 084

One IDEXX Drive • Westbrook, Maine 04092 USA

*Samples of this test kit model were independently evaluated by the AOAC Research Institute and were found to perform to the producer's specifications as stated in the test kit's descriptive insert. The producer certifies that this kit conforms in all respects to the specifications originally evaluated by the AOAC Research Institute as detailed in the PERFORMANCE TESTED certificate number 030302.

*SNAP and SNAPshot are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: idexx.com/patents.

© 2012 IDEXX Laboratories, Inc. All rights reserved. • 06-05061-03

IDEXX